

Logic[®] Vue Class II Enclosures

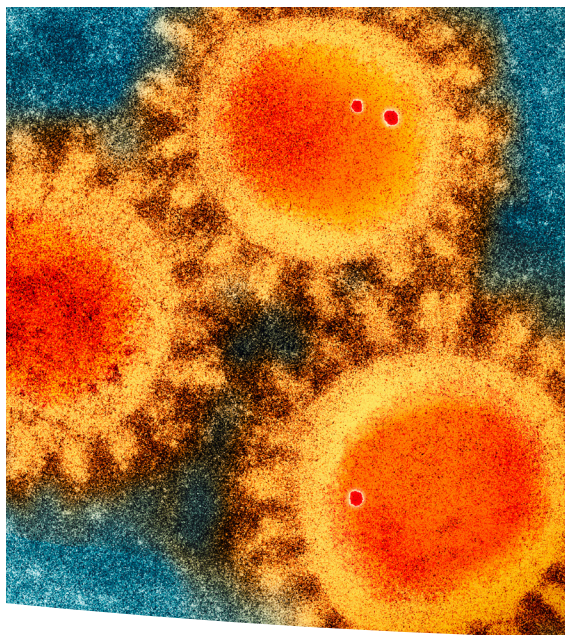




Table of Contents

Why the Logic Vue?	3	Ordering Information and Accessories	9
Enclosures At-a-Glance		Dimensions	10
Steps to Selecting a Biosafety Enclosure	4	Specifications	11
Biosafety Enclosure Product Lines	5		
Features and Benefits	6-8		

Go Big on biosafety.

Safety. Class II containment that meets NSF Standard 49 for personnel and product protection, plus robust construction ensure the Logic Vue keeps you safe while automating research.

Convenience. At 65% larger than a standard BSC, the Logic Vue is compatible with a variety of automated instruments. Accessibility and visibility from multiple sides allows for effortless operation and observation.


Intelligence. MyLogic™ OS is simple to operate and includes smart systems like, CAP™ technology, which continually monitors and auto-adjusts system airflows for maximum safety. Intuitive connectivity talks directly to automated instruments, pausing work when conditions are unsafe.

Don't trade safety for convenience. Choose both.



Labconco Biosafety Cabinets & Enclosures

Steps to Selecting a Biosafety Enclosure

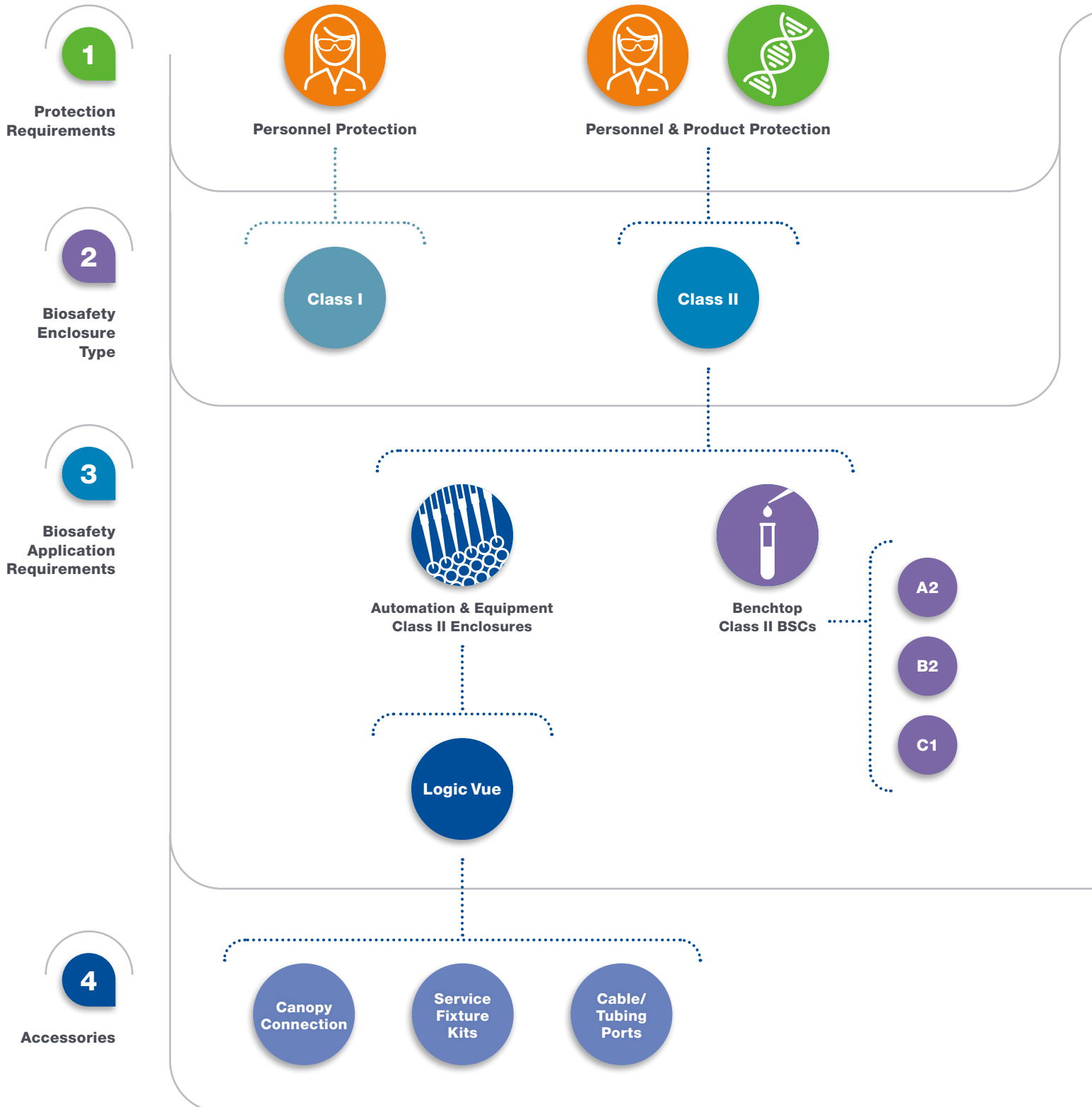

Labconco Biosafety Cabinets & Enclosures

Logic® and Axiom® Class II Products

Large Class II Biosafety Enclosures



Logic Vue



Class II Biosafety Cabinets



Logic+ Type A2




Logic+ Type B2



Logic Vue Accessories



Service Fixture Kits

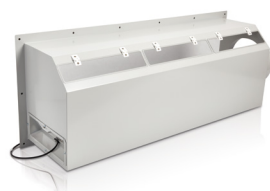

Canopy Connection



Cable/Tubing Ports




Axiom Type C1



Logic® Vue Class II Enclosures

Features

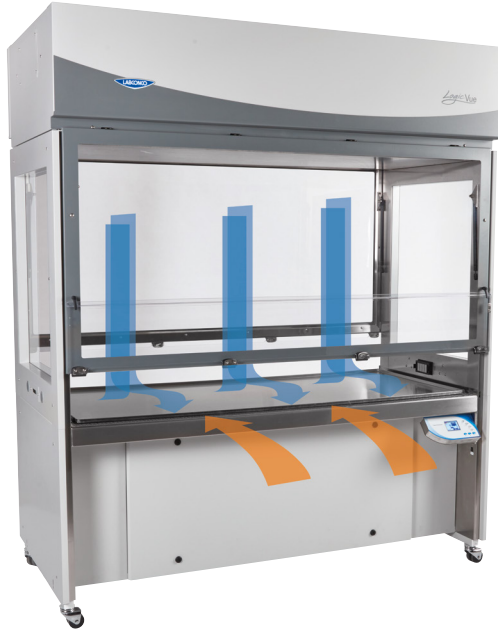

The Logic Vue Class II Enclosure

provides personnel and product protection:

- Meets NSF/ANSI Standard 49 for personnel and product protection
- Laminar ISO 5 or better air (99.99% efficient at 0.3 microns)
- Constant Airflow Profile™ auto-adjusting blowers contain biohazards while maintaining aseptic conditions within the enclosure




Convenience for Instrumentation

- Excellent visibility from all sides of enclosure
- Large 42.2" (107.2 cm) interior height provides ample space for tall instruments
- Deep 34.7" (88.1 cm) interior
- Spacious 31" (78.8 cm) interior working depth provides clearance for a variety of instruments
- Dual full-open sashes provide easy interior access. Work from the front and load or service equipment from either side.

Logic® Vue Class II Enclosures


Features

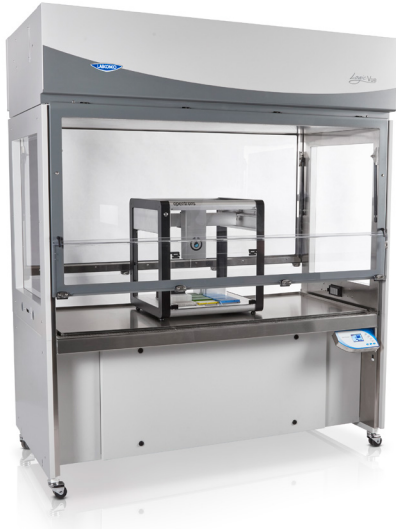

Convenient design provides unrestricted front-side access to entire work area during operation. An ergonomic knee well provides comfort while seated.



MyLogic™ OS intuitively communicates important status updates, such as alerts and filter life remaining. An intelligent safety interlock pauses automated instruments when operating conditions are unsafe.

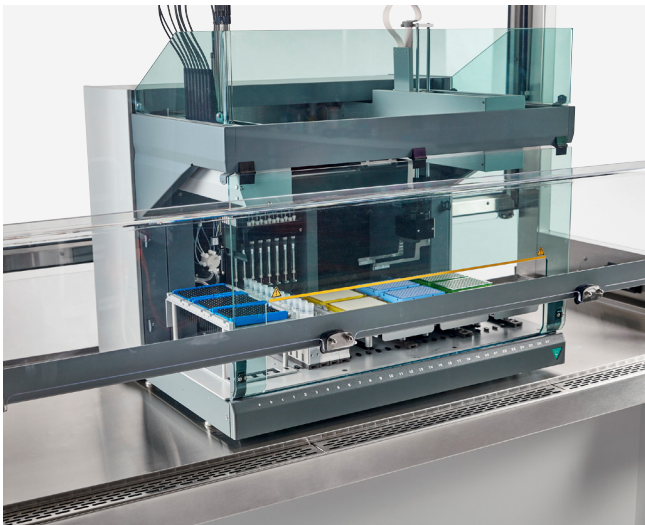

Robust work surface with included SpillStopper tray supports equipment up to 250.0 lbs (113.4 kg) per every two feet (61.0 cm).



Multiple pass-throughs allow safe passage of cables, cords, service fixtures, aerosol management systems, chillers and more.

Logic® Vue Class II Enclosures

Features & Benefits



Designed for instrumentation

including liquid handlers, cell sorters, and more.

Meets NSF/ANSI 49

requirements for personnel and product protection.

Ergonomic design

Ergonomic design allows setup for standing or sitting positions with built-in knee space.

HEPA supply and exhaust filters are minimum 99.99% efficient (at 0.3 microns), providing ISO 5 or better air.

CAP™ Technology continuously monitors and adjusts airflow to keep users safe.

Dual full-open sashes provide easy access for equipment loading and servicing.

Front sash opening provides unobstructed access to entire work surface during operation.

Stainless steel interior

Type 304 construction with reinforced work surface for easy decontamination.

Versatile design

Maximum visibility, pass-throughs, access doors, and monitor & keyboard arm compatible.

Ships fully assembled

Simplified installation with modular design to fit through standard doorways.

MyLogic™ OS Information Center



LCD Display

Clearly visualizes cabinet conditions in eight languages.

Filter Life Indicator

Accurately displays remaining HEPA filter life. Alerts users at 20% and 10% before alarming at 0%.

Blower Status

Displays when blowers are turned on, in reduced flow mode, or off.

Alerts & Alarms

Visual and audible signals notify users of unsafe cabinet conditions.

Safety Interlock

Optional connection pauses/stops automation equipment when sash is opened.

Data Connectivity

Exports system information via USB or dry contact relays for remote monitoring.

Logic® Vue Class II Enclosures

Ordering Information & Accessories

Catalog Number Configurator Use this key to configure the seven digit catalog number.

3	3	0	0	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
Select the WIDTH of your enclosure. This is the fifth digit of your catalog number.				Select the ELECTRICAL CONFIGURATION of your enclosure. This is the seventh digit of your catalog number.		
4 = 4' 6 = 6'				0 = 115V, 50/60 Hz, 30A		
Select the SHIPPING CONFIGURATION of your enclosure. This is the sixth digit of your catalog number.				1 = 230V, 50/60 Hz, 20A		
0 = Standard 1 = International Packaging*						

All models require (not included):

*Requires High Cube container.


- Dedicated electrical circuit

Shipping Data

Model	Shipping Width	Shipping Depth	Shipping Height	Shipping Weight
4'	67.5" (171.5 cm)	51.0" (129.5 cm)	104.0" (264.2 cm)	956 lbs (434 kg)
6'	91.5" (232.4 cm)	51.0" (129.5 cm)	104.0" (264.2 cm)	1275 lbs (578 kg)

Note: International Packaging height is 101.0" (256.5 cm)

Accessories

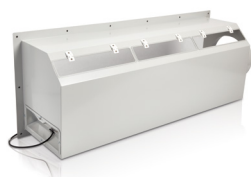

For thimble exhausting the Logic Vue. Canopy maintains a negative pressure connection to house exhaust. Detects insufficient exhaust in less than one second and signals MyLogic OS to display an alarm. Adds 11.8" (30.0 cm) to overall depth. Powder-coated

steel. Collar is 12" (30.5 cm) ID. Exhaust connection is required. **Exhaust damper is recommended and remote exhaust is required (not included). Installation required.** Shipping weight 80 lbs. (36 kg)


3347811 — 4' Logic Vue Canopy Connection.

Requires up to 1452 CFM @ up to 0.35" w.g.

3347813 — 6' Logic Vue Canopy Connection.

Requires up to 1953 CFM @ up to 0.35" w.g.

3349700—Safety Interlock Cable. Provides interface between the Logic Vue and automation devices that employ a safety interlock. Opening the sash of the Logic Vue triggers the interlock, immediately signaling a device to pause or stop. Shipping weight 3.0 lbs. (1.4 kg)

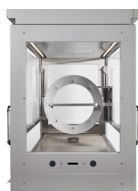

Large Access Port(s).

Provide extra access through sashes or side windows. Up to two per Logic Vue. Factory-install only. Contact Labconco for information.



3339000—Logic Vue Monitor & Keyboard Arm. Attaches to left side of Logic Vue. VESA mount supports monitors up to 18 lbs. **Installation is required.** Shipping weight 23 lbs. (10 kg)

3349800—Airflow Sensor Kit. Displays inflow and downflow velocities on MyLogic OS screen. Includes sensor, wiring harness and hardware. **Installation and certifier calibration required.** Shipping weight 3.0 lbs. (1.4 kg)



Large Tubing Pass-Throughs. For passage of larger tubing through sidewall of Logic Vue for external equipment such as aerosol management systems or chillers. Constructed of 16 gauge 304 stainless steel. Adds 1.8" (4.6 cm) to overall width. **Installation is required.** Shipping weight 3.0 lbs. (1.4 kg)

3349200—Large Tubing Pass-Through, 3/4" 0.75" OD (1.91 cm).

3349201—Large Tubing Pass-Through, 1-1/4" 1.25" OD (3.18 cm).


3768906—Replacement Prefilter. Extends life of the supply HEPA filter. Shipping weight 3.0 lbs. (1.4 kg)



3747500—Service Fixture Kit. Includes serrated hose tip with ball valve, quarter turn control handle and hardware. **Installation required.**

Logic® Vue Class II Enclosures

Dimensions

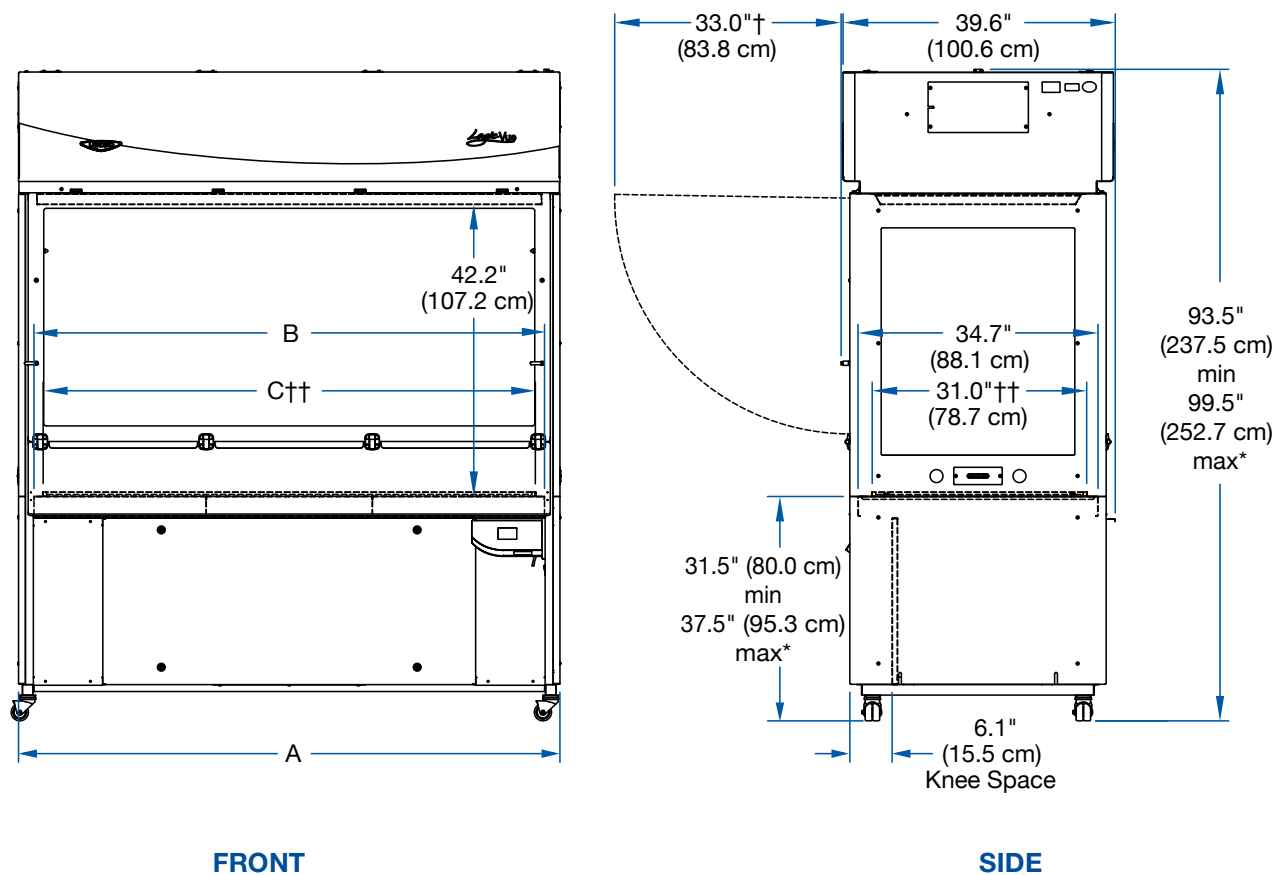

Model	Overall Width (A)	Maximum Interior Width (B)	Working Width (C)
4'	54.2" (137.7 cm)	49.8" (126.5 cm)	47.2" (119.9 cm)
6'	78.2" (198.6 cm)	73.8" (187.5 cm)	71.2" (180.8 cm)

*Adjustable in 1" (2.5 cm) increments.

†Swing clearance for Front Hatch. Same for Rear Hatch.

††Footing of instrumentation cannot exceed these dimensions. Maximum interior dimensions are 34.7" (88.1 cm) depth x Working Width (C).

NOTE: Logic Vue may be reduced in size to 77.5" (196.9 cm) height x 35.0" (88.9 cm) depth for passage through small doorways. Some disassembly required.

Logic® Vue Class II Enclosures

Specifications

All models feature:

- Intrinsically safe, single-pass negative pressure design
- Fully-opening front/rear sashes and side windows constructed of durable polycarbonate
- * 8.0" (20.3 cm) front lower sash opening with unrestricted access to full working area
- Unidirectional ISO Class 5/Grade A air or better
- HEPA filters, supply and exhaust, are 99.99% efficient or better on particles 0.3 microns or larger
- Individual HEPA filters are 100% scannable
- Standard pre-filter extends life of supply HEPA filter
- Spacious working area of 42.2" h x 31.0" d (107.2 x 78.7 cm)
- Loading height of 42.2" (107.2 cm)
- * MyLogic™ OS displayed on a full-color LCD screen with adjustable angle and touchpad controls
- * Smart-Start™ package automates blower and light functions for start-up and shut-down
- Energy-efficient electronically communicated motors (ECM)
- * Constant Airflow Profile (CAP) safe airflow monitoring system
- Electronic security lock
- * Robust multi-piece work surface accommodates equipment up to 250.0 lbs. (113.4kg) per every 2 feet (60.1 cm)
- 304 stainless steel interior
- Chemically-resistant epoxy powder-coated steel exterior
- * SpillStopper work surface tray
- * Ergonomic deep knee well
- * Adjustable work surface height of 31.5-37.5" (80.0-95.3 cm) in 1" increments (2.5 cm)
- * Remote monitoring via standard USB and dry contact relays
- * LED lighting illuminates work surface for maximum visibility of 133 footcandles (1432 lux)
- * Safety Interlock port for automated equipment
- * 16 cable/cord pass-through ports standard
- * 2 service fixture provisions per sidewall
- * 2 large pass-through provisions per sidewall
- * Ergotron® computer/media arm attachment point
- * Four transport casters and four seismic leveling feet
- Reduces to 77.5" h x 35.0" w (196.9 x 88.9 cm) for passage through small doorways
- * Ships fully assembled with minimal installation required
- One year warranty on parts and labor

* Labconco exclusive feature.

Standards conformance & regulations:

- Meets or exceeds NSF/ANSI Standard 49 for personnel and product protection
- CE Mark (230V models)
- UL Standard 61010-1
- CAN/CSA 22.2 No. 61010-1
- ADA-compliant operation

Options include:

- 115V or 230V electrical configurations
- International packaging configuration
- Factory-installed large access port(s)

All models require (not included):

- Dedicated hard-wired electrical circuit (select based on model)
 - 100-115V, 50/60Hz: 30A circuit
 - 208-230V, 50/60Hz: 20A circuit

Optional accessories:

- Canopy Connection Kit
- Monitor and Keyboard Arm
- Large Tubing Pass-Throughs (up to 2 per side)
- Service Fixtures (up to 2 per side)
- Airflow Sensor
- Safety Interlock Cable

**Visit labconco.com to request information
for all Labconco products.**

LABCONCO CORPORATION

8811 Prospect Avenue
Kansas City, MO 64132

(800) 821-5525 | (816) 333-8811

labconco.com



©2023 LABCONCO CORPORATION

Printed in the U.S.A.

Product design subject to change without notice.

2-32-2/15/23

